Soil Sampling Protocol eDentity [ENG]

Date: 27 June 2025

Purpose: Sampling soil for molecular analysis by the eDentity facility

Required Tools (Bring Your Own)

- Smartphone with camera and internet access
- Small shovel or spade (as clean as possible)

Required Tools (Provided in this sampling kit)

- 9 pre-labeled 15 mL sample tubes for DNA analysis
- Transect rope (5 m, with a 'knot' marking the 2.5m middle)
- Metal sampling spoon

Procedure

1. Select a Sampling Site

The site should be homogeneous in land cover and representative for the general area.

2. Mark the Sampling Area

Place the transect rope (5m) on the ground. This will mark the top of the transect location (sites 1, 2, 3 in figure 1a). Take out three random tubes and put them on the ground at the beginning, the middle (knot) and end of the transect line. Remove the transect rope and place it at a 90° angle from the tube at site 1. This will mark the sampling locations for points 1, 4 and 7 (figure 1a). Take out two tubes and put them on the ground at the middle and end of the transect line. Repeat this action to determine the location for sampling points 5 and 6 and 8 and 9 (figure 1a). If the area does not allow for a 5 m x 5 m grid, collect as many samples as possible without resizing the grid (figure 1b). In some cases, this may mean collecting only one sample. Always position the grid to ensure sample 1 can be collected (see figure 1b).

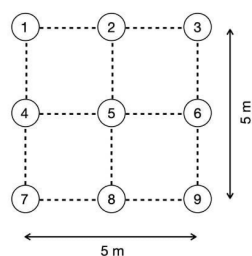
Figure 1a
sampling grid

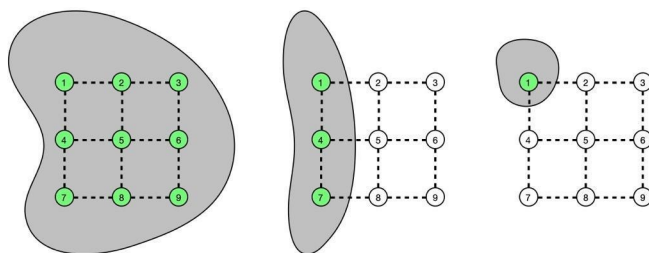
Figure 1b
sampling grid examples with insufficient area for sampling

3. Record Sampling Information

1. Open your browser and go to <https://ada.arise-biodiversity.nl/> or scan the QR code on the sampling kit.
2. Login with your username and password.
3. Click on the “+” and “eDentity” to enter sample information.
4. You now see the sampling grid on your screen. Click on the first sample in the upper left corner. Scan the barcode of the corresponding tube with your camera. If the barcode cannot be scanned, enter the barcode number manually.
5. Repeat the step above for all additional sampling points (if any). If all samples are added, press “next”.
6. Add a project number, location name (optional), and collector name. Add any labels that are relevant for this sampling plot (optional). Press “next”.
7. Location data (date, time, position) are auto filled. Make changes if necessary. Hit “next”.
8. Select the relevant land use and soil type categories from the dropdown menu. Add kitnumber (ED-SOIL-XXXX) and other remarks if necessary. Hit “next”.
9. Take a picture of the sampling grid showing all grid corners, and with sampling point 1 in the upper left corner of the picture. Press “next”.
10. You now see an overview of the sampling data. Press “Save” when no changes need to be made. You are now ready to collect the samples.

use your Google
Chrome browser

4. Collect Soil Samples

1. Take **sample tube 1** from sampling point 1.
2. Remove any litter (loose material) from sampling point 1.
3. Take a spade or small shovel and put it in the ground (~10 cm) at point 1.
4. Bring the soil to the surface and fill the tube **no further than the mark** (red line) with multiple small scoops of soil. Avoid large roots or stones as much as possible.
5. Repeat this step for all remaining sampling points. Cleaning the spade or shovel between sampling points is not required.

5. Avoid Cross-Contamination

- If sampling a new location, use a fresh eDentity sampling kit.
- Do not reuse any items from the kit.
- Clean the spade or shovel as much as possible with tap water.

6. Storage and shipping instructions

- Send the sample kit including ALL contents back within 3 days after sampling using the supplied Reply Number.
- If shipping within 3 days is not possible, store the kit in the refrigerator at 2-8°C for up to 14 days.

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